

The Design and Development of an Automated Glycemic Index Analyser

Phillip Clancy, Next Instruments Pty Ltd, Condell Park, NSW, Australia.



Glycemic Index: Health Issues

NutriScan

Glycemic Index of the foods we eat is important for the following reasons:

- 1. Diebetes the control of blood glucose and Insulin
- 2. Obesity appetite control, appetite suppression
- 3. Children energy and obesity
- 4. Heart Disease weight control
- 5. Colon Rectal & Bowel Cancer- efficient gastric flow





Glycemic Index Analysis


- NutriScan
- •Glycemic Index measures the release of glucose from food under enzymatic digestion.
- •GI relates mostly to foods that contain measureable levels of carbohydrates: cereals, dairy, baked goods, energy bars, etc
- Glucose has a GI of 100
- GI ranges from:
 - High GI: > 70
 - Medium GI: 55-69
 - Low GI: < 55



Glycemic Index Analysis



- Gold Standard In Vivo Method
- -10 subjects -12 hour fast, no exercise
- Eat 50 g of food sample
- Blood glucose measured each 15mins for 2 hours
- 3 tests: 1 glucose reference

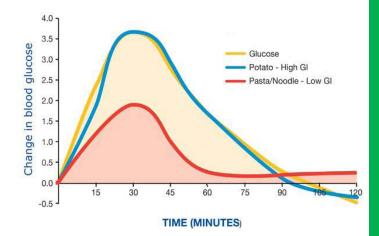

2 repeat tests for food samples

- Area under the curve is the Glycemic Response GR
- GI = 100* (GR for 50g of food/GR for 50g of Glucose)



NutriScan GI and RS Analysers



Developed by CSIRO, Div of Food and Nutrition Sciences

- GI = Glycemic Index
- RS = Resistant Starch
- 1st in the world
- Artificial Gut to simulate the digestion of food and to measure the glucose released over:

5hr (GI Analysis) and 16hr(RS Analysis) period.






NutriScan GI and RS Analysers

- The system involves a series of incubations, at physiological pH and temperature that essentially mimic the buccal, gastric and pancreatic phases of digestion.
- Intended for use in Research into Animal and Human Nutritional.
- Intended not as Replacement for In Vivo method.
- Intended for Product Development
 and QC





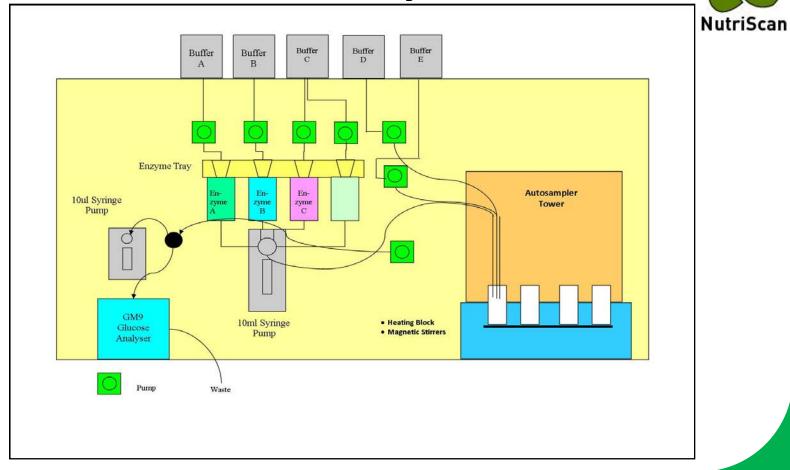





Pre-weighed Enzyme Cartridge

- 5, 10 and 20 sample cartridges
- Ensures consistency of results against the validation studies performed by CSIRO
- Reduces errors due to weighing







Pumping System

- 4 peristaltic pumps
 - 1 x 3 head pump loads enzyme bottles
 - 2 single heads pumps load other buffers
 - 1 single withdraws samples from digest.
- 2 syringe pumps
- 10ml syringe dispenses enzyme solutions
- 500ul syringe injects 10ul sample into GM9





NutriScan

20 Cup Digester

- XYZ Rotary Dispensing System
- Fully programmable
- Gentle Agitation :
 20 position magnetic
 stirrer block
- Temperature controlled heating block: 37C





Glucose Analyser

Glucose Oxidase Reaction

- Measures the depletion of Oxygen using an Oxygen Electrode
- RS232 coms to external PC








PC Controlled

- Fully programmable
- Program
 - Enzyme sequencing
 - Incubation times
 - Sampling times
 - Buffer loading volumes and sequences.
- Automatic data report
 - Plots Glycemic Load,Hydrolysis Index
 - Reports GI



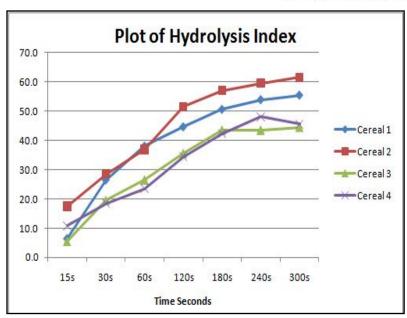



Comparison between the In Vivo GI and RS VS the In Vitro methods:

NutriScan	


Parameter	In Vivo Method GI RS		In Vitro Methods GI RS	
Time required To produce results	7 days	24 days	1 day	1 day
Cost per test	\$4000	\$8000	\$25-100	\$25-100
Sample throughput (Runs per week)	2	1-2	100-200	50
Discrimatory Power	15GI units	0.5g RS/g food	5 GI units	0.2g RS/g food
Inter Assay Precision CV ²	30-40%	20-45%	2-10%	5-20%



CSIRO Food and Nutrition Sciences have analysed hundreds of foods using their GI and RS Analysers and compared the results with the In Vivo Methods.



Sample Type	GIValue	CV%
Shortbread Biscuit	69	3.4
Graham Cracker	83	2.7
Savoury Cracker	80	3.8
Breakfast Cereal 1	82	4.9
Breakfast Cereal 2	52	2.3
Breakfast Cereal 3	65	2.7
Snack Food	75	4.4
Fruit Bar	44	10.9
White Bread	82	3.8
Wholemeal Bread	74	2.7

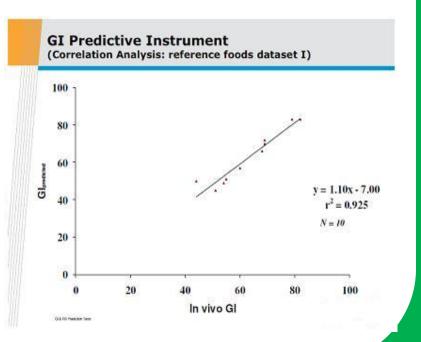




Resistant Starch is the starch that passes through to the small intestines without undergoing enzymatic digestion.



Sample Type	RSValues
Pasta Breakfast Cereals Bread - Wholegrain Bread - White Potato - Fries - Powdered Pasteries Legumes	.5-1.0% .2-3.0% 1.5-2.5% .5-1.0% 1.0-1.5% .3% .39% 2.0-5.0%

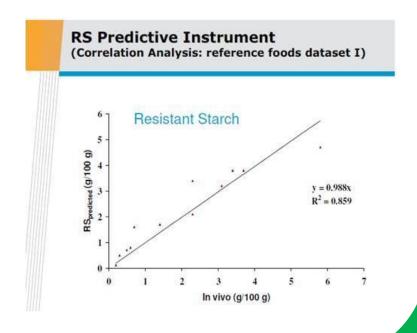



Summary:



NutriScan GI20 Glycemic Index Analyser

- Fully Automated Artificial Gut
- 20 Sample Capacity
- Improved Precision vs In Vivo Method
- Validated by CSIRO
- Suitable for Product Development and Routine QC

Next Instruments, Booth No. 7355

